



Assembly of RBR S14


For assistance, please contact info@spinchem.com.

SpinChem AB
Tvistevägen 48C
90736 Umeå
Sweden


Page 1. Loading the RBR S14
Page 16. Disassembly for cleaning
Page 19. Reassembly for storage

1. Loading the RBR S14

Loading the RBR S14

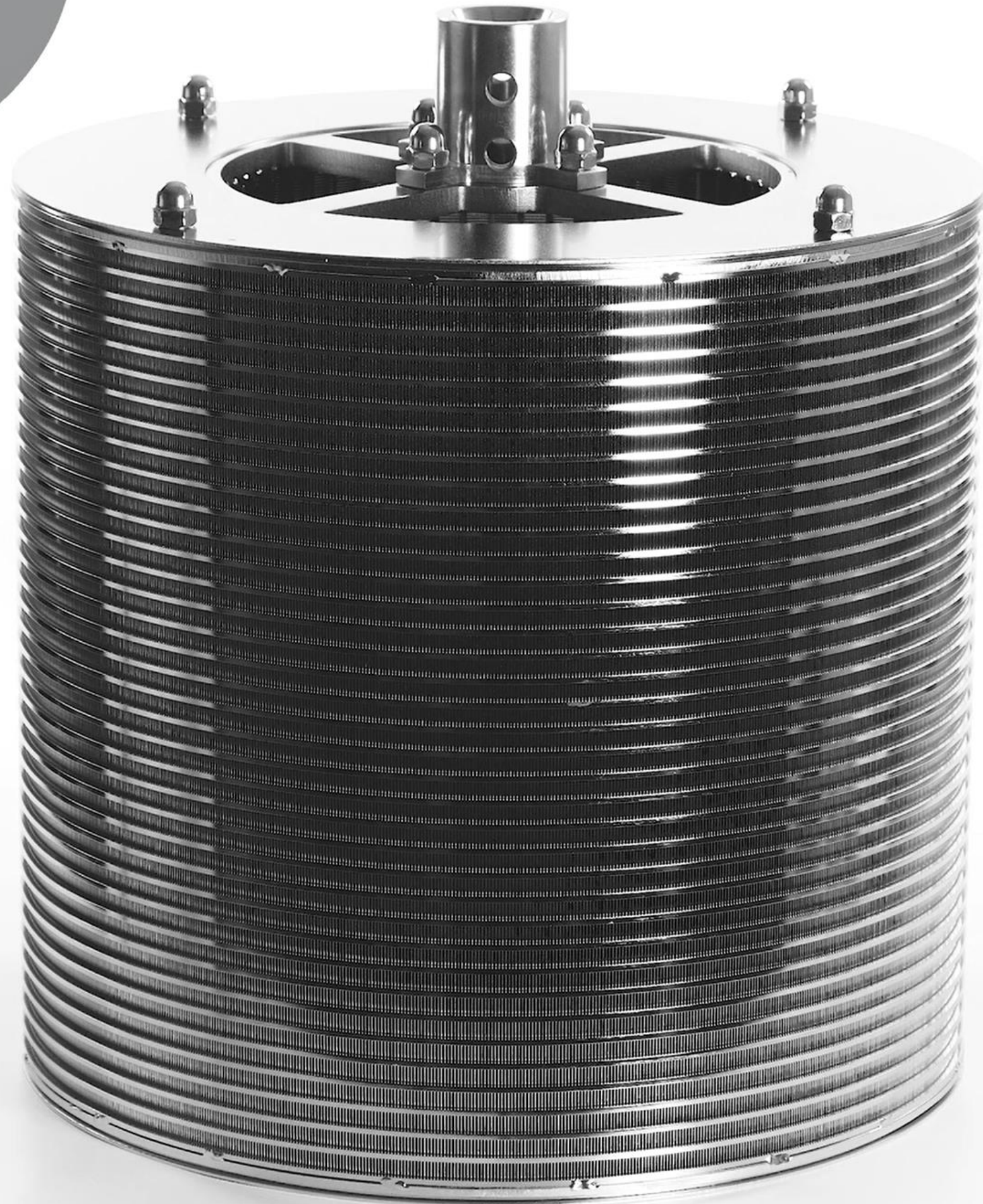


Loading the RBR S14

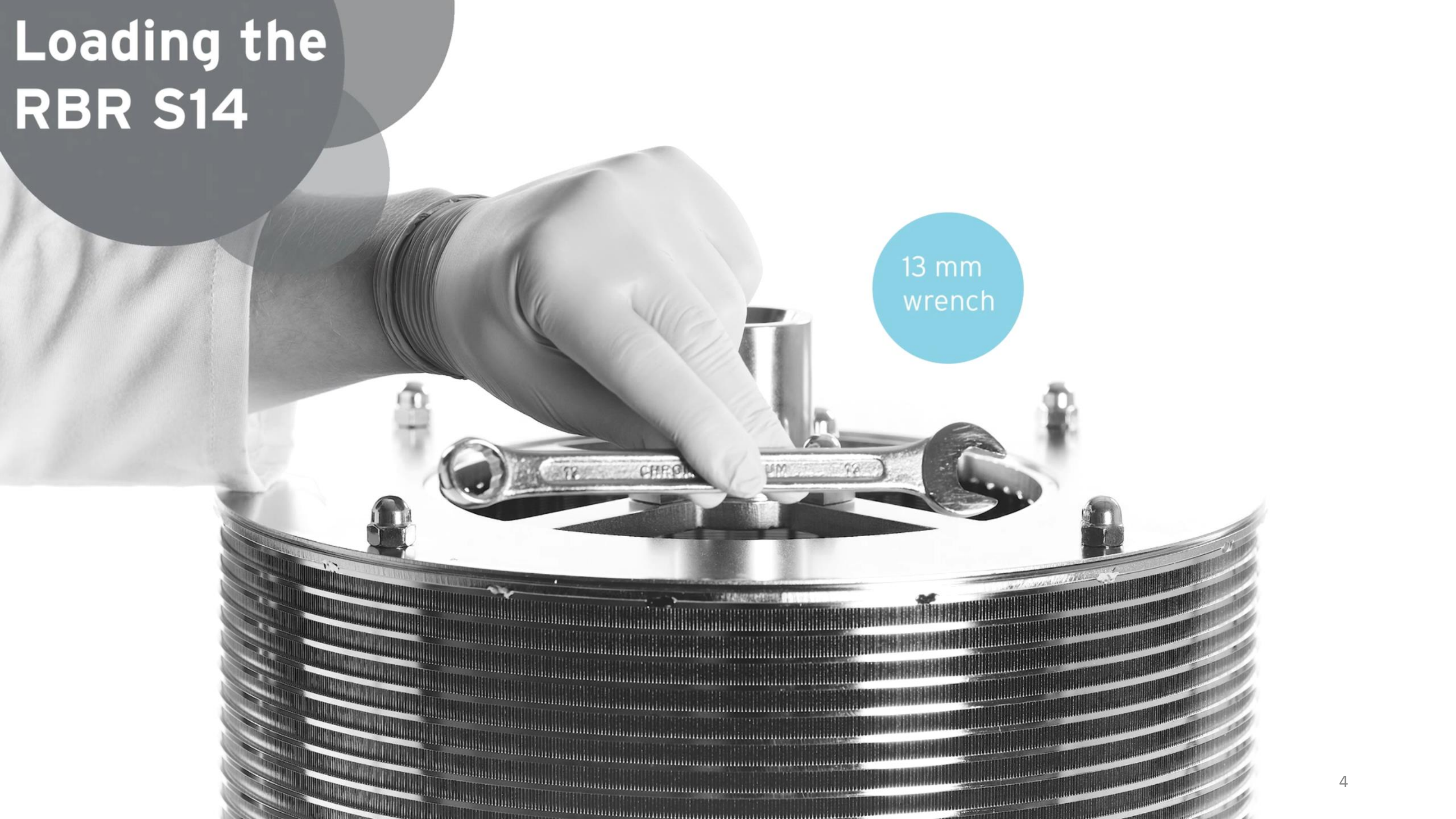

13 mm
wrench

Loading the RBR S14

Release the four nuts on the lid of the RBR S14

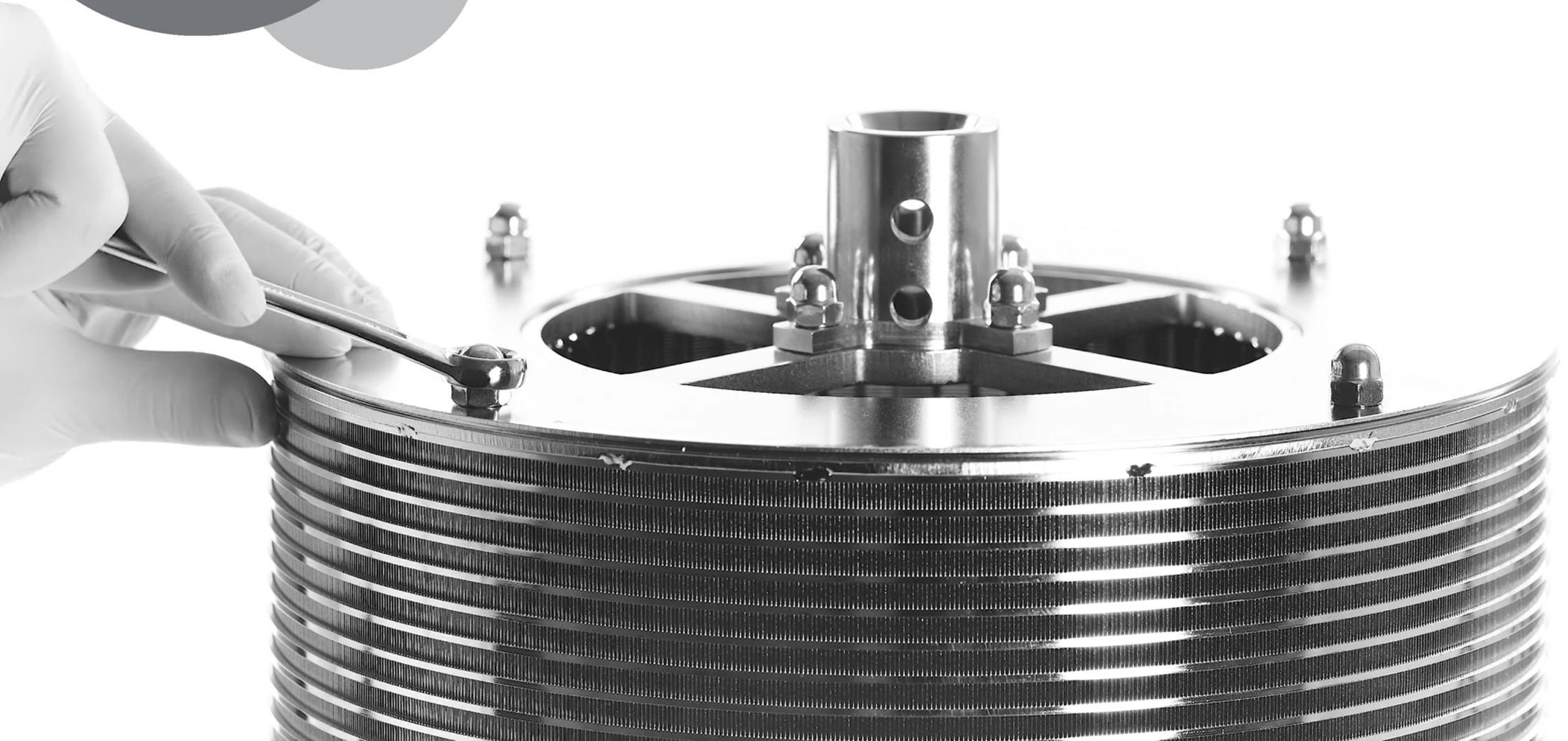

Loading the RBR S14

Pry the lid open with your
hands or a plastic tool



Loading the RBR S14

Remove the lid



Loading the RBR S14

1st option:
solid phase
free in RBR



Fill the RBR with free solid phase directly from a container, bag, etc.

Loading the RBR S14

2nd option:
solid phase
baged in RBR




Fill the RBR with bags (sold separately by SpinChem AB), packed with solid phase.

Loading the RBR S14




Loading the RBR S14

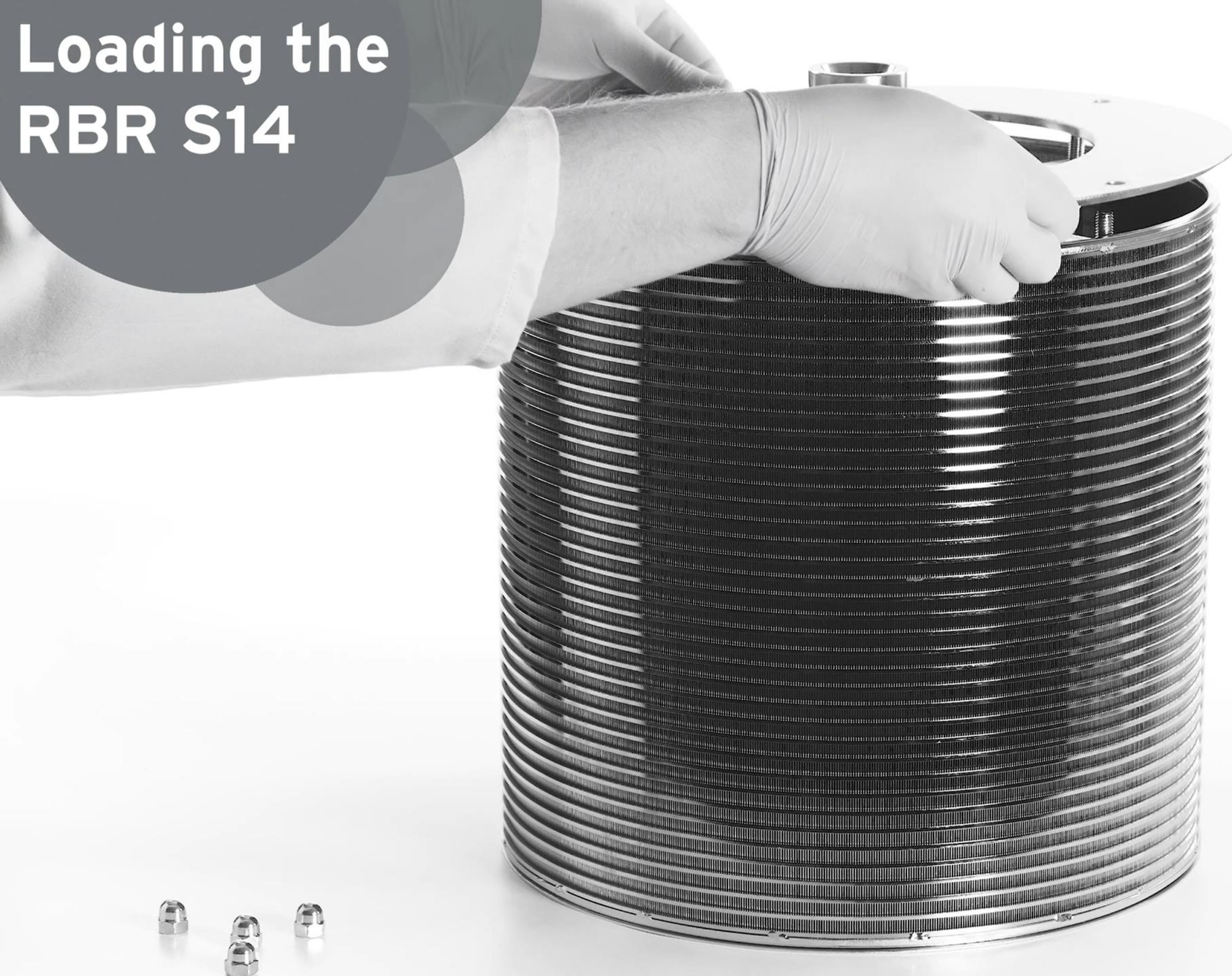

Align the lid with the
threaded rods on the RBR
S14

Loading the RBR S14


Firmly press down on the lid,
ensuring that it seals against the
outer and inner filter cylinders

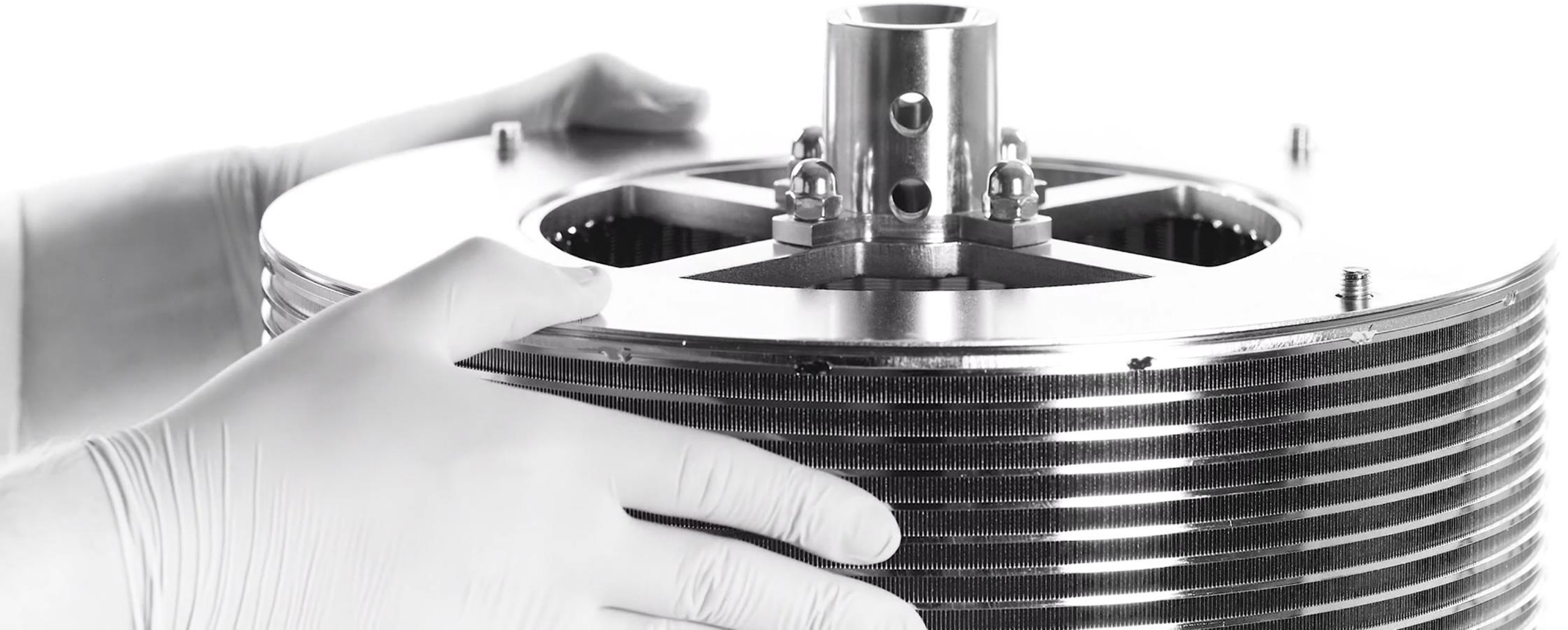

Loading the RBR S14

ensure that the lid is seated all the way

Firmly press down on the lid, ensuring that it seals against the outer and inner filter cylinders

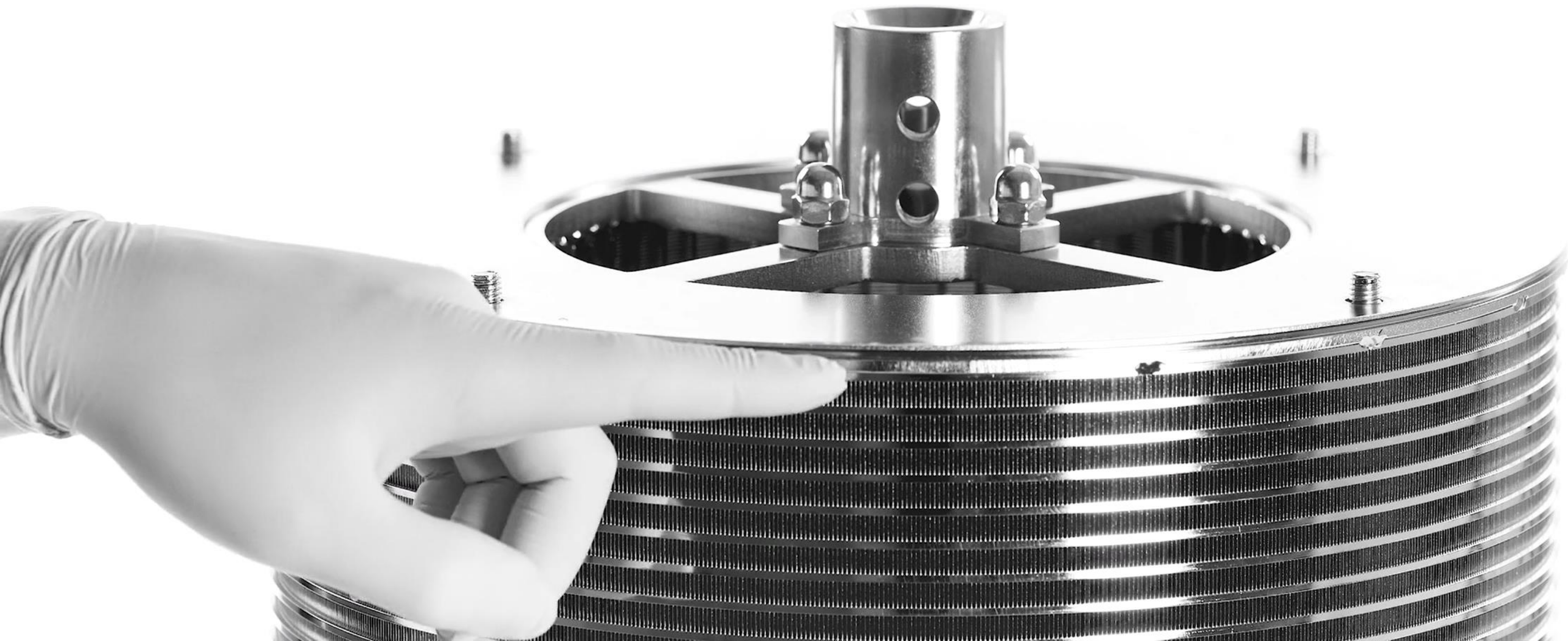

Loading the RBR S14


ensure that the lid is seated all the way

Note: Align the number on the lid with the corresponding number on the RBR



Loading the RBR S14

Fasten the four nuts on the lid of the RBR S14

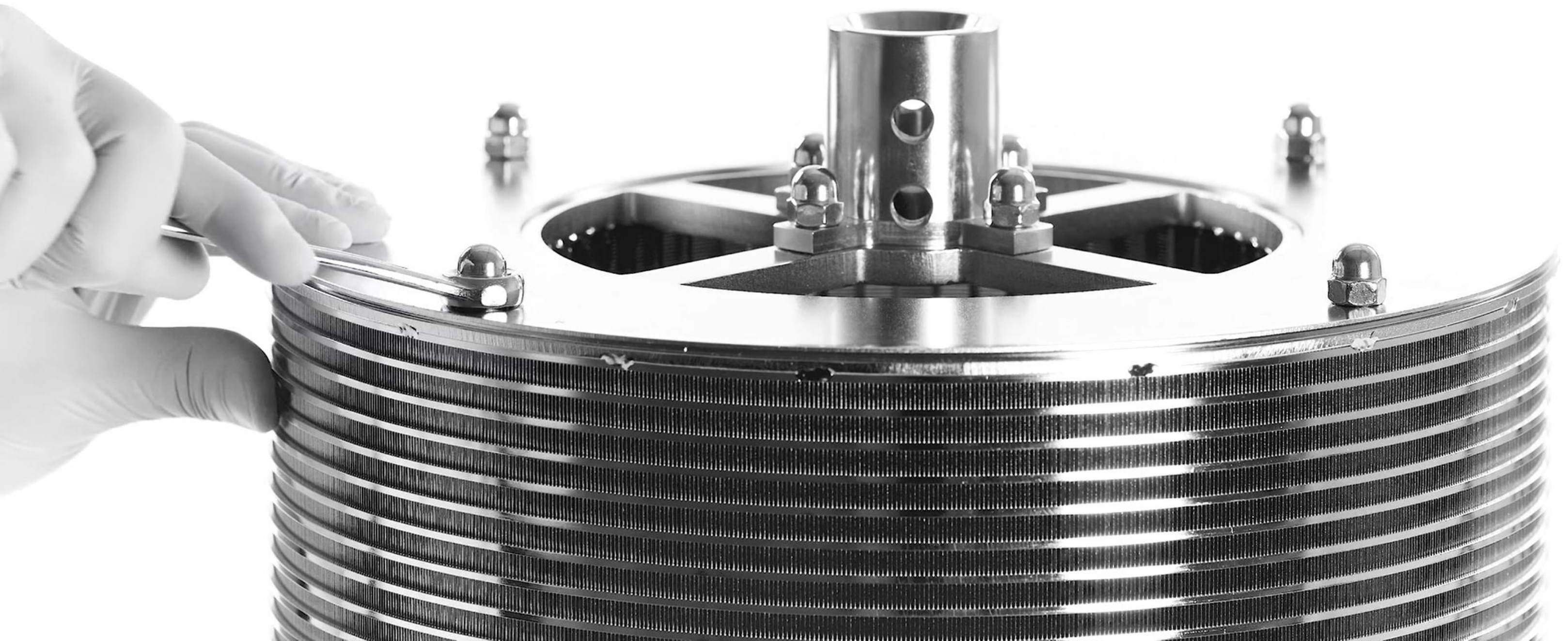

2. Disassembly for cleaning

Disassembly for cleaning



Grip the outer filter cylinder with your fingers and press down on the inner cylinder with your thumbs

Disassembly for cleaning

Cleaning of the filter cylinders is most efficiently done using pressurized air, water or solvent jets or with a plastic bristled brush



3. Resassembly for storage

Reassembly for storage



Place the outer filter
cylinder over the inner
filter cylinder

Reassembly for storage



Firmly press down on the outer filter cylinder, ensuring that it seals against the bottom plate

Reassembly for storage

ensure that
the filter
is seated all
the way



Firmly press down on the
outer filter cylinder,
ensuring that it seals
against the bottom plate

Reassembly for storage



Align the lid with the
threaded rods on the RBR
S14

Reassembly for storage




ensure that
the lid is
seated all
the way

Firmly press down on the lid, ensuring that it seals against the outer and inner filter cylinders

Reassembly for storage

ensure that
the lid is
seated all
the way

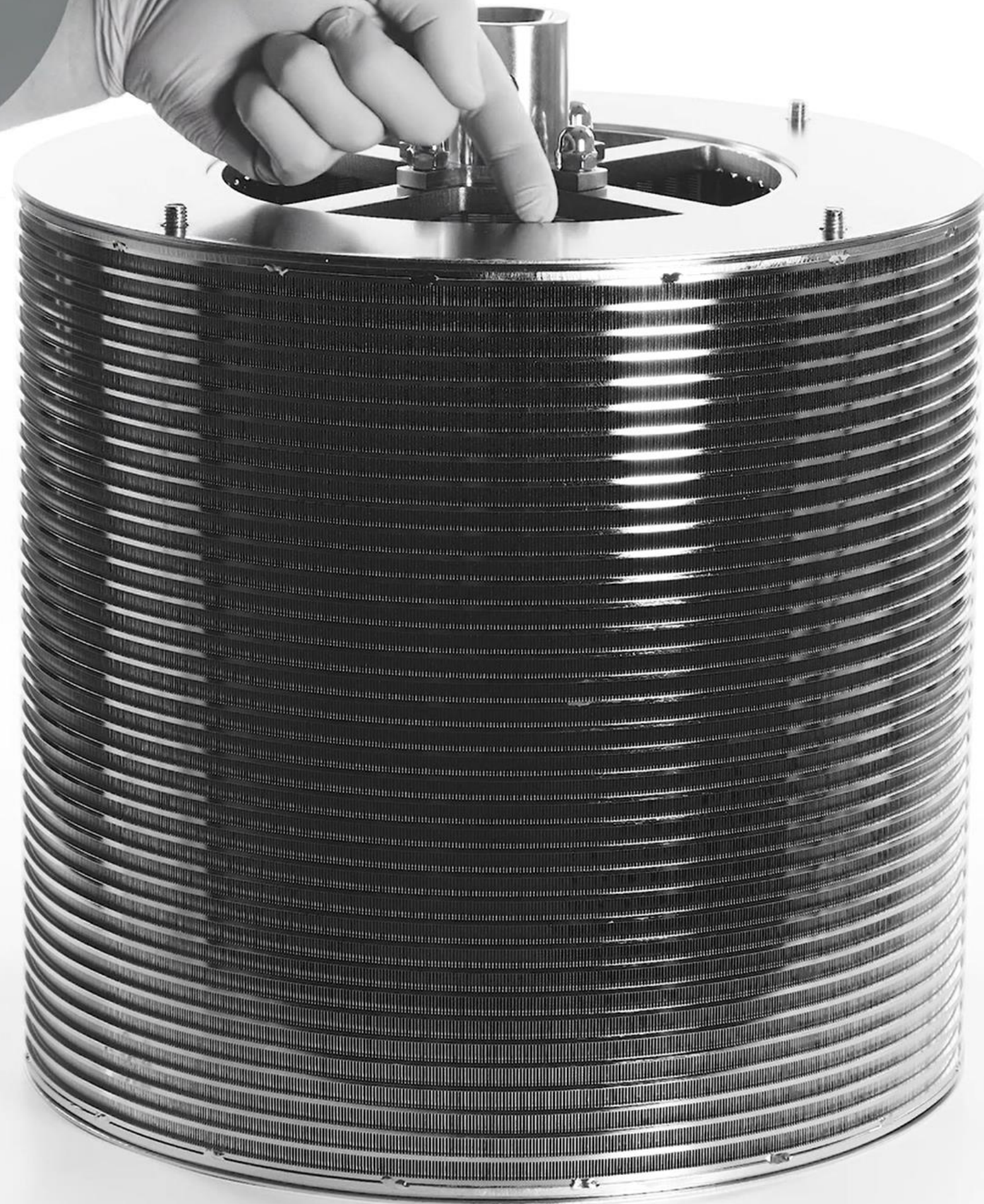

Firmly press down on the lid, ensuring that it seals against the outer and inner filter cylinders

Reassembly for storage

Fasten the four nuts on the lid of the RBR S14

